



PCR BASED QUARANTINE PROGRAM

(e.g. one line of n=10 Transgenic mice) Each mouse is sampled directly.

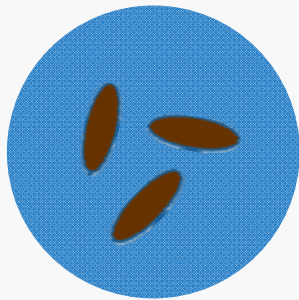


Fecal pellet

Collect one fecal pellet from each mouse, pool up to 10:1 into a single tube

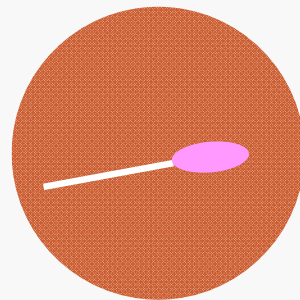


Body swab

Use a single sticky swab to cover the entire pelt of the animal, including perianal hairs, pool swab tips up to 10:1 into a single tube

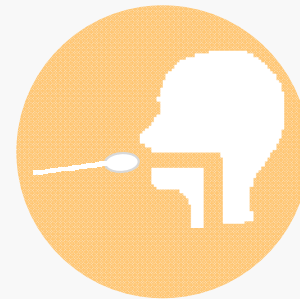

Oral swab

Use a small cotton swab to sample the mouth of each mouse, pool swab tips up to 10:1 into a single tube

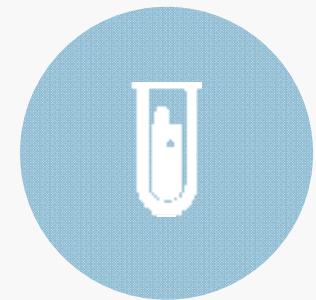

One PRIA PCR sample

Samples will be pooled into a single test article for testing

* Include survival serology to determine previous exposure